

**CAMPYLOBACTER SELECTIVE SUPPLEMENT, CCDA**

FOR 500 ml OF CAMPYLOBACTER BLOOD-FREE (CCDA) AGAR

Freeze-dried mixture for the isolation of *Campylobacter* spp.

Description	
Code number:	4 ml: CCS80004
Colour:	White
Appearance:	Homogeneous lyophilisate

**Direction:** Suspend **24 g of Campylobacter Blood-Free (CCDA) Agar Base (CCA20500)** in 500 ml of distilled water and heat with frequent agitation until the medium boils well. Sterilise by autoclaving at 121 °C for 15 minutes. Cool to 50 °C and add aseptically the contents of **one vial of Campylobacter Selective Supplement, CCDA (CCS80004)** reconstituted with 4 ml of sterile distilled water. Mix well before pouring.

**FORMULA [mg/vial]**

Cefoperazone	16
Amphotericin B	5

**Storage conditions:** Protected from light, at 2-8 °C.

**Warning!**

Use before the expiry date on the label.

**In vitro diagnostic – for professional use only!**