

COLIFORM SELECTIVE SUPPLEMENT

FOR 500 ml OF ChromoBio® COLIFORM ES

Freeze-dried mixture for the simultaneous detection of coliforms and *Escherichia coli*.

Description	
Code number:	4 ml: COS80004
Colour:	White
Appearance:	Homogeneous lyophilisate

Direction: Suspend **15 g of ChromoBio® Coliform** in 500 ml of distilled water and heat with frequent agitation until the medium boils well. Sterilise at 100 °C (in water bath or flowing steam) for 30 minutes. Cool to 50 °C and add aseptically the contents of **one vial of Coliform Selective Supplement (COS80004)** reconstituted with 4 ml of sterile distilled water. Mix well before pouring.

FORMULA [mg/vial]

Cefsulodin	2,5
Vancomycin	2,5

Storage conditions: Protected from light, at 2-8 °C.

Warning!

Use before the expiry date on the label.

In vitro diagnostic – for professional use only!