

LISTERIA SELECTIVE SUPPLEMENT, OXFORD

FOR 500 ml OF LISTERIA SELECTIVE AGAR, OXFORD

Freeze-dried mixture for the isolation of *Listeria monocytogenes* according to ISO 11290-1.

Description	
Code number:	4 ml: LS080004
Colour:	White
Appearance:	Homogeneous lyophilisate

Direction: Suspend **29,5 g of Listeria Selective Agar Base, Oxford (LA020500)** in 500 ml of distilled water and heat with frequent agitation until the medium boils well. Sterilise by autoclaving at 121 °C for 15 minutes. Cool to 50 °C and add aseptically the contents of **one vial of Listeria Selective Supplement, Oxford (LS080004)** reconstituted with 4 ml of 1:1 mixture of ethanol and sterile distilled water. Mix well before pouring.

FORMULA [mg/vial]

Cycloheximide	200,0
Colistin	10,0
Fosfomicin	5,0
Acriflavine	2,5
Cefotetan	1,0

Storage conditions: Protected from light, at 2-8 °C.

Warning!

Use before the expiry date on the label.

In vitro diagnostic – for professional use only!