

TTC SOLUTION, STERILE

FOR THE PREPARATION OF 10 OR 30 LITRE OF MEDIA

TTC (2,3,5-Triphenyl-2*H*-tetrazolium chloride) solution for the preparation of some media.

Description	
Code number:	10 ml: TTC80010, 30 ml: TTC80030
Colour:	Water clear
Appearance:	Transparent solution

Direction: Prepare 500 ml of sterile medium base from one of the dehydrated medium bases for the detection of enterococci or coliforms (e.g. CATC Agar Base, KF Streptococcus Agar Base, Slanetz-Bartley Agar Base, Tergitol 7 Agar Base) according to the direction of the given medium. Cool to 50 °C and add aseptically **10 drops (0,5 ml) of TTC Solution, Sterile (TTC80030)**. Mix well before pouring.

FORMULA

TTC	5 %
Solvent	Distilled water

Storage conditions: Protected from light, at 2-8 °C.

Warning!

Use before the expiry date on the label.

In vitro diagnostic – for professional use only!