

TRIBUTYRIN SUPPLEMENT

FOR 500 g OF TRIBURYRIN AGAR BASE

Tributyryn (1,2,3-Tributyrylglycerol) for preparation of Tributyrin Agar.

Description	
Code number:	250 ml: TRS80250
Colour:	Water clear
Appearance:	Transparent

Direction: Suspend **20 g of Tributyrin Agar Base (TRA20500)** in one litre of distilled water. Add **10 ml of Tributyrin Supplement (TRS80250)** and mix until homogeneous. Heat with frequent agitation until the medium boils well. Sterilise by autoclaving at 121 °C for 15 minutes. Cool to 50 °C with frequent agitation and pour plates immediately to solidify quickly.

Warning!

The ready medium must be homogeneous turbid gel!

FORMULA

Tributyryn (bacteriological grade)	250 ml
------------------------------------	--------

Storage conditions: Protected from light, at room temperature.

Warning!

Use before the expiry date on the label.

In vitro diagnostic – for professional use only!