TECHNICAL SHEET



VIOLET RED BILE GLUCOSE AGAR, PH EUR

A glucose containing selective and differential medium for the detection and enumeration of Enterobacteriaceae according to PH EUR (Agar Medium F – Crystal Violet Neutral Red Bile Agar with Glucose).

Dehydrated media		
Code number: 500 g: VBE20500, 5 kg: VBE25000		
Colour:	Beige	
Appearance:	Homogeneous hygroscopic powder	
pH before sterilization (25 °C):	7,2 – 7,6	

Direction: Suspend **51,5 g** in one litre of distilled water and heat with frequent agitation until the medium boils well. Mix well before pouring.

Warning!

The medium is heat sensitive. No further sterilisation is necessary or desirable.

Prepared media			
Bottled media:	100 ml: VBE30100, 500 ml: VBE30500		
Plated media:	55 mm: VBE50055, 90 mm: VBE50090		
Colour:	Reddish purple		
рН (25 °C):	7,3 – 7,5		

Direction: Dispense the melted bottled media aseptically into sterile Petri-dishes. Media in Petri-dishes are ready to use.

FORMULA in g/l

Gelatin peptone	7,000
Yeast extract	3,000
Bile salts	1,500
Glucose monohydrate	10,000
Lactose monohydrate	10,000
Sodium chloride	5,000
Neutral red	0,030
Crystal violet	0,002
Agar	15,000

Note: The typical formula can be adjusted to obtain optimal performance.

Storage conditions: Store the dehydrated media tightly closed in a dry place at room temperature. Store the bottled media protected from light at room temperature. Store the plated media protected from light at 2-8 °C. Use before the expiry date on the label.

Quality control:

Test strains	Incubation temp: 37 °C	Growth	Incubation time: 24 h
Escherichia coli	ATCC 25922	Good, purplish red colonies with precipitate halo	
Proteus mirabilis	ATCC 29906	Good, purplish red colonies without swarming	
Enterococcus faecalis	ATCC 29212	Inhibited	

References: European Pharmacopoeia

In vitro diagnostic - for professional use only!